## INSPECTION REPORT

MODEL \_\_\_\_\_

SRT1

### ELECTRICAL SAFETY

1. Earth continuity
2. Insulation
3. Flash test

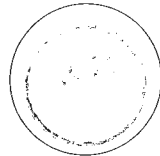
### FUNCTIONAL

1. Indicators
2. Roller control
3. Visual acceptance

QUALITY CONTROL INSPECTOR

## Bibby Sterilin

### Bibby Sterilin Ltd

Beacon Road Stone  
Staffordshire ST15 0SA  
United Kingdom  
Tel: +44 (0)1785 812121  
Fax: +44 (0)1785 813748  
www.bibby-sterilin.com

### Bibby Sterilin France

BP 79-77793 Nemours Cedex  
France  
Tel: +33 1 64 45 13 13  
Fax: +33 1 64 45 13 00  
bibby@bibby-sterilin.fr

### Bibby Sterilin Italia

Via Alcide de Gasperi 56  
20077 Riozzo Di Cerro Al Lambro  
Milano  
Italy  
Tel: +392 983 8969  
Fax: +392 982 30211

### Afora S.A.

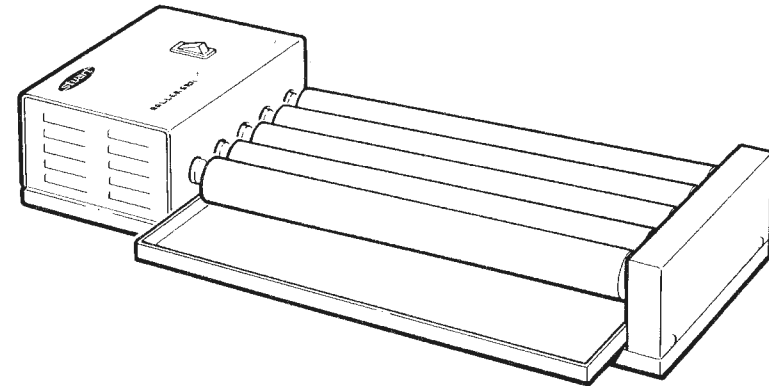
Calle Aribau 740  
08006 Barcelona  
Spain  
Tel: +343 93-306 98 000  
Fax: +343 93-306 98 2 3  
e-mail: marketing@aloua.com  
www.afora.com

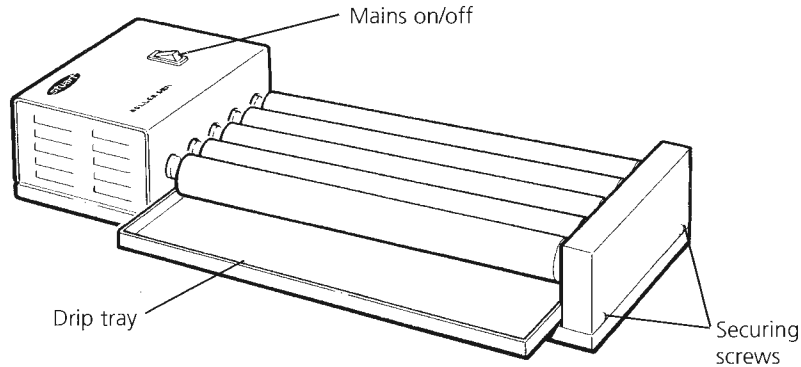
## Bibby Sterilin

Roller Mixers  
SRT1 & SRT2

Instructions for use  
Manuel d'utilisation  
Istruzioni per l'uso  
Instrucciones de funcionamiento  
Bedienungsanweisung

**Thank you for purchasing this piece of Bibby Sterilin equipment. To get the best performance from the equipment, and for your own safety, please read these instructions carefully before use.**

This equipment is designed to operate under the following conditions:-

- ❖ For indoor use only
- ❖ Use in a well ventilated area
- ❖ Ambient temperature range +5°C to +40°C
- ❖ Altitude to 2000m
- ❖ Relative humidity not exceeding 80%
- ❖ Mains supply fluctuation not exceeding 10%
- ❖ Overvoltage category II IEC60364-4-443
- ❖ Pollution degree 2 IEC664

If the equipment is not used in the manner described in this manual the protection provided by the equipment may be impaired.

## Electrical Installation

**THIS EQUIPMENT MUST BE EARTHED**

**Before connection please ensure that the line supply corresponds to that shown on the rating plate. This model requires a supply rated at 220 - 240V, 50 Hz, ~, single phase.**

### Power requirements

Model	Wattage
SRT1& 2	50W

There is an IEC socket at the rear of the instrument for connection to the mains supply.

The unit is supplied with two mains leads fitted with IEC plugs for connection to the instrument.

One has a UK 3 pin plug and the other has a 2 pin "Shuko" plug for connection to the mains.

Choose the lead appropriate for your electrical installation and discard the other. Should neither lead be suitable, take the lead with the UK plug and replace the plug with a suitable alternative.

See the enclosed instruction sheet for advice on how to carry out this procedure.

Should the mains lead require replacement a cable of 1mm<sup>2</sup> of harmonised code H05W-F connected to an IEC 320 plug should be used.

### IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN

The mains lead should be connected to the instrument BEFORE connection to the mains supply.

## General Description

The Stuart Roller Mixers provide a gentle, but highly efficient, rocking and rolling action. The rollers rotate at a fixed speed of 33rpm and rise and fall 16mm.

Most sizes of tube, Bijoux, Universals and other bottles can be accommodated. For large capacity bottles (up to 120mm diameter), one or more of the rollers is easily removed, using the spanner provided, to give improved grip. Similarly, it is possible to quickly convert the unit to provide a 'roll only' action to function as a simple roller.

**SRT1:** Has 5 rollers

**SRT2:** Has 10 rollers

## Safety Advice Before Use

- ❖ The unit should be carried using both hands with fingers under each side frame.
- ❖ Never move or carry the unit when in use or connected to the mains electricity supply.
- ❖ Do not mix flammable liquids or use the equipment in hazardous atmospheres.
- ❖ In case of mains interrupting, a fault or mechanical failure, the unit will continue to operate on removal of fault.
- ❖ Mechanical energy can lead to breakage of glass vessels. Use with care.

## Operation

Position the apparatus on a firm level surface and connect to the mains electricity supply. Place the samples to be mixed between the rollers. Once all the samples have been set, switch the Roller Mixer ON using the Mains On/Off switch.

**WARNING: Ensure the unit is switched Off at the mains before altering roller configuration.**

### Removing the rollers

The rollers may be removed from your unit to accommodate larger vessels. To remove, first loosen the two securing screws shown in figure 1 and remove the end panel.

Undo the pivot bolt of each roller to be removed using the spanner provided. Lift the free end of the roller and gently pull it away from the drive unit. Keep all rollers and bolts in a safe place.

### Roll only operation

Your SRT Roller Mixer is factory set for a 'rock and roll' mixing motion. The action can be converted to a roll only operation.

First, remove all rollers as described above. One by one, starting at one end, replace each roller using the **CENTRIC** securing hole in both the drive and end pivots. Tighten the bolts using the spanner provided.

## Maintenance & Servicing

**WARNING: Ensure the unit is disconnected from the mains electricity supply before attempting maintenance or servicing.**

Periodically clean the instrument using a damp cloth and mild detergent solution. Do not use harsh or abrasive cleaning agents.

**Any repairs or replacement of parts MUST be undertaken by suitably qualified personnel.**

The following spares and accessories are available from your laboratory dealer.

Description	Catalogue Number
Roller	DSRT1014
Spanner	USRT1C2

For a comprehensive list of parts required by service engineers conducting internal repairs and a service manual, please contact the Technical Service Department of Bibby Sterilin Ltd. quoting both the model and serial number.

Only spare parts supplied or specified by Bibby Sterilin Ltd. or its agents should be used. Fitting of non-approved parts may affect the performance and safety features designed into the instrument.

If in any doubt, please contact the Technical Service Department of Bibby Sterilin Ltd. or the point of sale.

Bibby Sterilin Ltd  
Stone  
Staffordshire ST15 0SA  
United Kingdom  
Tel: +44 (0) 1785 812121  
Fax: +44 (0) 1785 813748  
E-mail equipmnet@bibby-sterilin.com  
www.bibby-sterilin.com

## Technical Specification

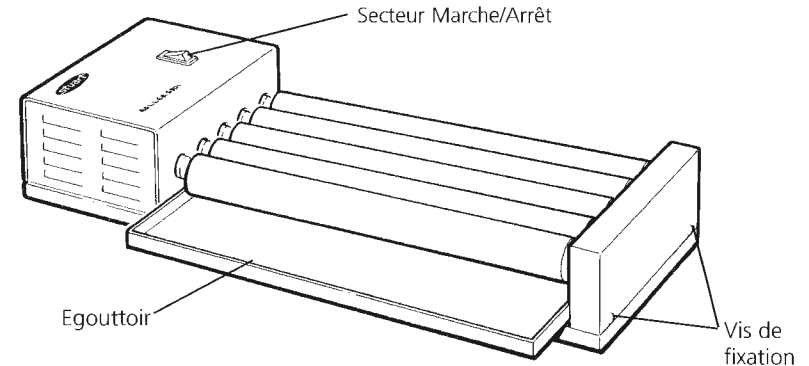
	SRT1	SRT2
No. of rollers	5	10
Speed	33rpm	33rpm
Amplitude	16mm	16mm
Roller size (length x diameter)	340 x 30mm	340 x 30mm
Overall dimensions (WxDxH)	500 x 200 x 90mm	500 x 380 x 90mm
Net weight	5kg	7kg

## Warranty

Bibby Sterilin Ltd warrants this instrument to be free from defects in material and workmanship, when used under normal laboratory conditions, for a period of two (2) years. In the event of a justified claim Bibby Sterilin will replace any defective component or replace the unit free of charge.

This warranty does NOT apply if damage is caused by fire, accident, misuse, neglect, incorrect adjustment or repair, damage caused by incorrect installation, adaptation, modification, fitting of non approved parts or repair by unauthorised personnel.

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## Notes

These products meet the relevant EC harmonised standards for radio frequency interference and may be expected not to interfere with, or be affected by, other equipment with similar qualifications. We cannot be sure that other equipment used in their vicinity will meet these standards

and we cannot guarantee that interference will not occur in practice. Where there is a possibility that injury, damage or loss might occur if equipment malfunctions due to radio frequency interference, or for general advice before use, please contact the Technical Service Department of Bibby Sterilin Ltd.

### *Declaration of Conformity*

#### *Model SRT1 & SRT2 Roller Units*

These Stuart Scientific products comply with the requirements of the EU Directives listed below:

**89/336/EEC** Electromagnetic Compatibility Directive, amended by 92/31/EEC and 93/68/EEC.

**73/23/EEC** Low Voltage Directive amended by 93/68/EEC

Compliance with the requirements of these Directives is claimed by meeting the following standards:

**EN55014-1:1997** Limits and methods of measurement of radio disturbance, characteristics of electrical motor-operated and thermal appliances for household and similar purposes, electric tools and similar apparatus.

**EN55104:1995** Electromagnetic Compatibility - General Immunity Part 1. Residential, commercial and light industry.

**EN61010-1:1993** Safety requirements for electrical equipment for measurement, control and laboratory use.

The technical documentation required to demonstrate that the products meet the requirements of these Standards is available for inspection on request.

Signed:  (Mr C. R. Baggaley)

Date: 2/6/02

Authority: Technical Manager

for

Bibby Sterilin Ltd - Stone - Staffs - ST15 OSA - UK  
Tel: +44 (0) 1785 812121 - Fax +44 (0) 1785 813748